



# **AIVION StreamWindow**

## **USER MANUAL**

AIVION, Jahnstraße 12, 85661 Forstinning

Tel: +49(0)8121-2208-0 Fax: +49(0)8121-2208-22

[www.aivion.com](http://www.aivion.com) [sales@avion.com](mailto:sales@avion.com)

AIVION is a Brand from eVision Systems GmbH

This application note is a technical support document and not an official part of the product documentation. In case of deviations from the datasheet please use the datasheet as reference information. Make sure that your board revision fits to the referred revision in the datasheet.

# Table of contents

- 0. Additional information..... 3
- 1. Setup the application ..... 3
- 2. Menubar ..... 3
  - 2.1 Camera..... 3
  - 2.2 Devices..... 3
- 3. Statusbar ..... 4


## 0. Additional information

The AIVION StreamWindow software only works together with the USB interfaces (TL6032, TL6035, TL6636, USB3 Neo). This is because the video data must be transferred via a USB cable.

## 1. Setup the application

The live image from the camera is displayed automatically as soon as the correct camera is selected (see chapter 2.2).

## 2. Menubar

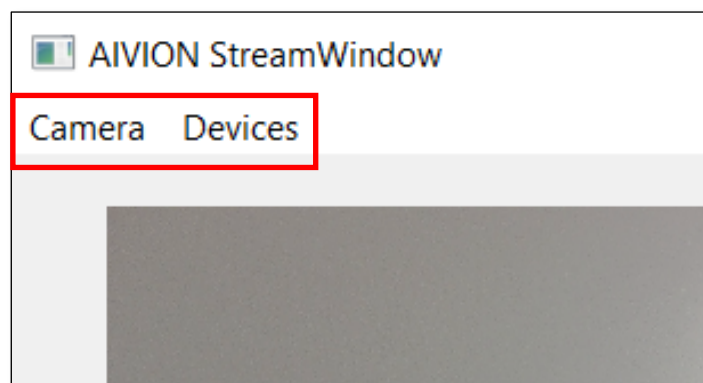

### 2.1 Camera

Under 'Camera' you can stop and restart the live image. With 'Capture Photo' you can take a still image of the live image. 'Exit' closes the application.

### 2.2 Devices

Under 'Devices' you can select the required camera from the connected cameras.

### 3. Statusbar



When you take a still image, the location where the image is saved is displayed.  
The default storage path for the screenshots is C:\Users\*your user*\Pictures.